



Pro-X Versaflo PPE Equipment

Instructions on how to put on Versaflo Safety Equipment and operate it.

Written By: Ami Manish Patel



Step 1 — Attaching the Filter



- Remove clear plastic lining from filter and line up filter with the outside of the Versaflo waist band
- Click in left side and then push in right side until it clicks.

Step 2 — Attach Filter Cover




- Place the light grey cover on top of the filter with the cut out towards the top (closest to hose hole) until you hear a click and all the sides are flush with the waistband.

Step 3 — Inserting Air Hose




- Place the hose that is attached to the head cover into the hole at the top of the waistband.
- Push down and twist clockwise to lock it into place.
- Double check it's locked by slightly tugging on it.

Step 4 — Adding Headpiece to Head Cover

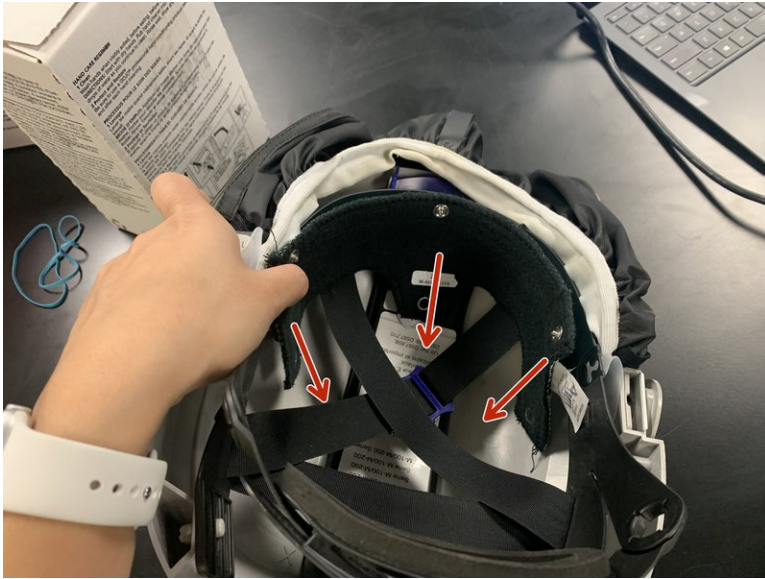

- Place the cloth headpiece with the blue side touching the top of the head cover.
- The three tabs should be facing down into the head cover.
- Wrap the three tabs under the plastic rim and button all three sides.

Step 5 — Attaching the Battery

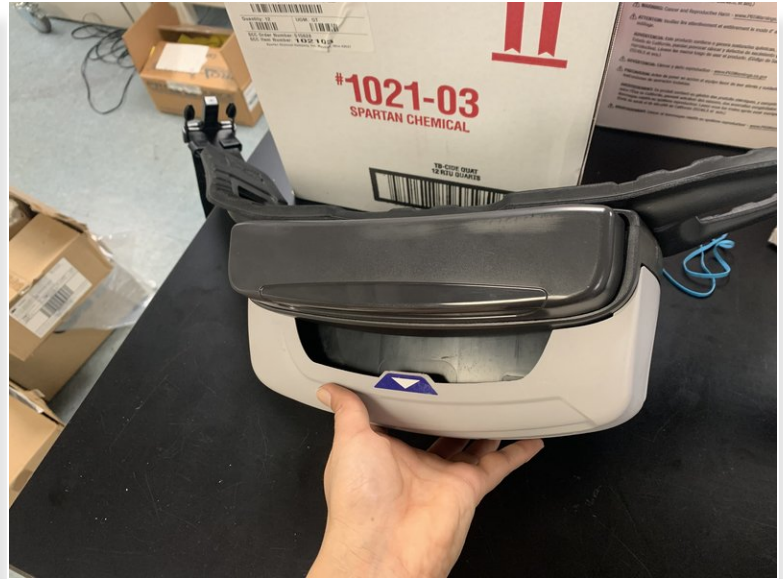

- Aligning the strip on the battery to the strip on the waistband
- Attach the battery pack to the bottom of the waistband until you hear the click

Step 6 — Starting the Air Flow



- Adjust the sizing on the head cover by twisting the knob and pulling the strap on the waistband. Strap them on.
- Hold down the power button on the top left corner for three seconds to start the air flow.
- Hold down power button for 3 seconds to turn off air flow once you are finished


This document was generated on 2022-01-01 01:57:45 AM (MST).

Step 7 — Adjusting Air Flow





- As you can see, the battery level and air flow levels are displayed at the top in green.
- In order to increase or decrease airflow, hold down the fan button that is next to the power button on top of the waistband.

Step 8 — Clean up

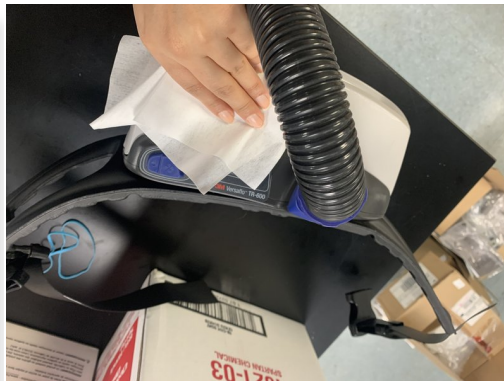

- When done using, remove headpiece, filter, cover, battery and hose.
- Use spray cleaner to wipe down headpiece and all surfaces before putting equipment back in cabinet.

Step 9 — Charge Battery

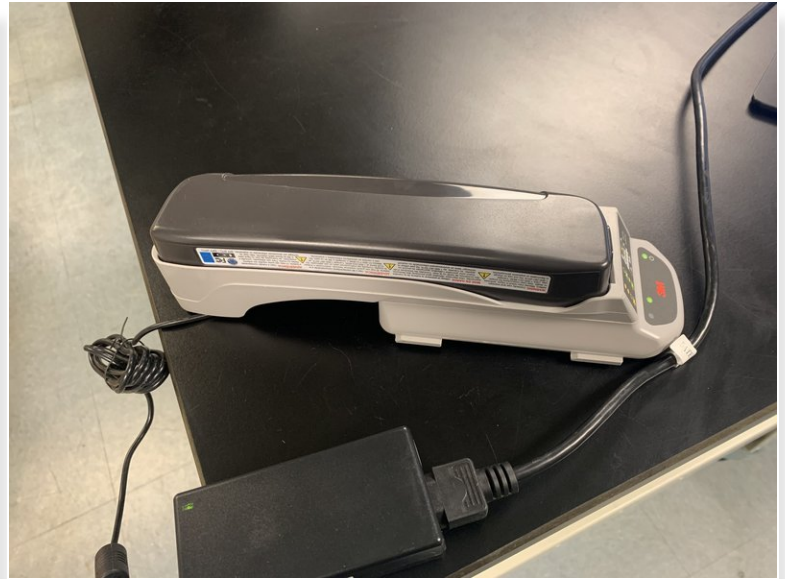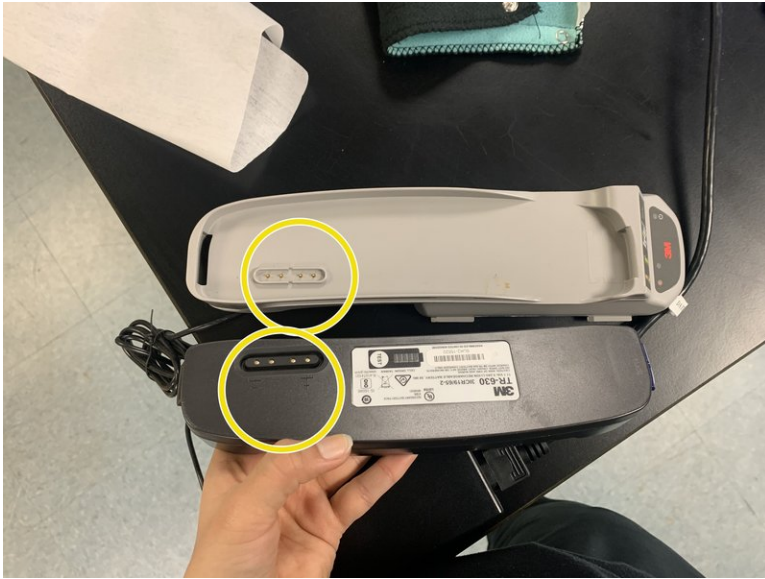

- Plug in battery charger to an outlet
- Match up battery and click into charger until all sides are flat and green light is on