

Keyence VR-3200 Executing a Basic 3D Scan

How to execute a basic scan using the Keyence

Written By: Amera Grace Gianotti



Step 1 — Running the Scan



- The magnification can be adjusted by selecting between low mag and high mag
 - Low mag is capable of scanning up to 1 cm depth, allowing for 12x, 25x, 38x, and 50x magnification
 - High mag is capable of scanning up to 1 mm depth, allowing for 40x, 80x, 120x, and 160x magnification
- The camera brightness can be adjusted, but **auto** is sufficient for most cases
- The focus can be adjusted manually using focus adjustment, or auto focus can be run with a double left click on the section you want to be focused
- After adjusting magnification and focus, run the scan by clicking measure

Step 2 — Stitching Multiple Scans



- You can only scan what is shown on the screen, so larger magnifications may require stitching in order for the object to be fully captured
 - Switch the stitching option to auto
- Under Area specification method, select Detailed. Underneath that, select Start point and end point
- Set the scanning boundaries by selecting four different positions of the object and assigning them to either top, left, right, or bottom
 - Use the arrows or click and drag to move the object on the screen to get the desired positions
- The boundaries you select are the outermost places the scan will be captured
- Click **Measure** when done setting the boundaries so the software runs the scan
 - The Viewer will automatically run the Analyzer software after the scan in complete

Step 3 — Exporting the Scan Result

- The scan result can be exported as a 3D CAD model
- Go to file, export, output 3D-CAD data
- Select stl file (.stl) as the file output file format